


Recommended procedure to mount syringe heaters:

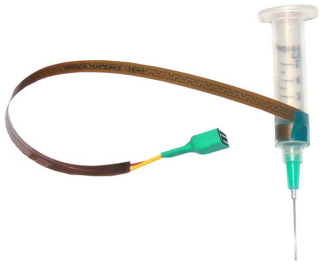

1. Position the heater on the syringe, the sensor down between the heater and the syringe surface, and fix the end in place using a small piece of an electrical tape.

2. Wrap the heater around the syringe, and fix the other end of the heater with the second piece of the electrical tape.



3. Make a loop using provided Velcro tape, and tighten it around the heater so the connector is sticking outside from the tape. Wrap the remaining Velcro tape around the syringe to prevent heat dissipation. You can use two Velcro tapes on both ends of the heater.

4. The surface of the larger volume syringes does not have to be tightly covered by the heater. Instead, the heater can spread along the whole surface.

